

Glacios

36k× Data Collection 1.125 Å/ubpix

LMM Settings

- Spot Size 6, 150 μm C2, LM 62×, C2 big enough to cover detector, Default Montage Settings

MMM Settings

- Spot Size 6, 50 μm C2, LM 380×, C2 70%

Low Dose Settings

- Spot size 6, 50μm C2
- View: NanoProbe, SA 6700×, 79% C2
- Record/T/F: NanoProbe, SA 36k×, C2 39.000%
- EPS 20.5, 1.125 Å/ubpix

Gain Ref Path

- C: drive>Program Data>Gatan>Reference Images
- File Format .dm4, most recent large file