

North Krios Tomography Mode 14eps, Pixel 3.44Å

LMM Settings

- Spot Size 7, C2 100µm, EFTEM LM 175×, Ill Area ~780µm, Default Montage Settings

MMM Settings

- Spot Size 7, C2 100µm, EFTEM Microprobe SA 3600×

HMM Settings

- Spot Size 7, C2 100µm, EFTEM Microprobe SA 3600×

Low Dose Settings

- Spot Size 7, C2 100µm
- View - EFTEM Microprobe SA 2250× - SA3600×
- Record - EFTEM Microprobe SA 26k×, Ill Area 3.3µm
- EPS ~14

Gain Ref Path

- C: ProgramData>Gatan>Reference Images>Most recent and Largest File size ~360,000KB
- File Format .DM4