

## North Krios Tomography Mode, 28eps, Pixel 3.44Å

### LMM Settings

- Spot Size 6, C2 100µm, EFTEM LM 175×, Default Montage Settings

### MMM Settings

- Spot Size 6, C2 100µm, EFTEM Microprobe SA 3600×, Ill Area ~40µm

### HMM Settings

- Spot Size 6, C2 100µm, EFTEM Microprobe SA 3600×

### Low Dose Settings

- Spot Size 6, C2 100µm
- View - EFTEM Microprobe SA 2250× - SA 3600×, Ill Area ~40µm
- Record - EFTEM Microprobe SA 26k×, Ill Area 3.3µm
- EPS ~28

### Gain Ref Path

- C: ProgramData>Gatan>Reference Images>Most recent and Largest File size ~360,000KB
- File Format .DM4